LymphoTranscript

ENGLISH	2
FRANCAIS	21

ENGLISH

GENEXPATH LymphoTranscript user guide

User precautions

RUO

This product is intended for molecular biology research purposes only. Do not use it for medical diagnosis, prevention or treatment purposes.

It is for professional use only.

Read all information in this user guide before use.

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contact@genexpath.com

support@genexpath.com

ENGLISH	2
Important precautions.....	5
General recommendations	5
Icons.....	5
Intended use	5
Test principle.....	6
Reagents	7
Content of GENEXPATH LymphoTranscript reagent kit.....	7
Format of reagent kits sold and quantities:.....	8
Reagents not supplied in the reagent kit:.....	8
Required equipment:	9
Before starting	9
Biological samples.....	9
Solutions to prepare	10
Programming the thermocyclers	10
Programme 1: Pre-PCR.....	10
Programme 2: PCR.	11
Detailed protocol	11
Step 1: Reverse transcription.....	11
Required reagents.....	11
Step 1.a: Reverse transcription; denaturation of RNA and hybridization of random hexamers.....	11
Step 1.b: Reverse transcription; conversion of RNA to cDNA.....	12
Step 2: Hybridization of probes	12
Required reagents.....	12
Hybridization of probes.....	13
Step 3: Ligation.....	13
Required reagents	13
Ligation	13
Step 4: Amplification and incorporation of barcodes and adapters	14

Required reagents	14
Amplification	15
Step 5: Purification and assay of sequencing libraries	15
Required reagents	15
Step 5.a: Purification of sequencing libraries	15
Step 5.b: Assay of sequencing libraries:.....	16
Step 6: Dilution, pooling and sequencing of libraries.....	16
Required reagents	16
Sequencing on an Illumina MiSeq sequencer	16
• Step 6.a: Dilution and pooling of libraries.....	16
• Step 6.b: Denaturation and dilution of the libraries pool.....	17
• Step 6.c: Preparation of the sequencing primer	17
• Step 6.d: Preparation of the injection sheet	17
• Step 6.e: Start of sequencing	17
Sequencing on an Illumina NextSeq 500/550 platform	18
• Step 6.a: Dilution and pooling of libraries.....	18
• Step 6.b: Denaturation and dilution of the libraries pool.....	18
• Step 6.c: Preparation of the sequencing primer	18
• Step 6.d: Preparation of the injection sheet	18
• Step 6.e: Start of sequencing	19
Step 7: Results analysis	19
Limits of the procedure	20
Bibliography	20
Symbol table	20
Notes.....	20

Important precautions

General recommendations

- For Research use only
- Follow laboratory best practices in terms of handling PCR products (wear disposable overalls and gloves, mark out dedicated zones for pre- and post-PCR, use filter tips).
- Also take precautions to avoid nuclease contamination likely to cause RNA and DNA degradation (use nuclease-free reagents and consumables).
- Ensure that the thermocyclers are in working order and calibrated based on manufacturer recommendations.
- It is particularly important not to substitute reagents not included in the kit, particularly buffers and enzymes used during the reverse transcription, ligation and PCR amplification steps. The incubation temperatures and times, as well as volumes and concentrations, must also be respected.
- **GENEXPATH LympHoTranscript** reagents are only intended for use with Illumina's MiSeq or NextSeq 500/550 sequencing platforms.
- Safety data sheets are available in the user space
- If the user lights out errors in the instruction manual: please send an email at contact@genexpath.com
- Any serious incident occurring in connection with the device must be notified to us at contact@genexpath.com.

Icons

Important points and critical steps of the protocol that could compromise result quality.

Steps where the protocol may be suspended.

Intended use

This protocol is intended for **GENEXPATH LympHoTranscript** testing. It is used to prepare sequencing libraries for Illumina's MiSeq or NextSeq 500/550 sequencers.

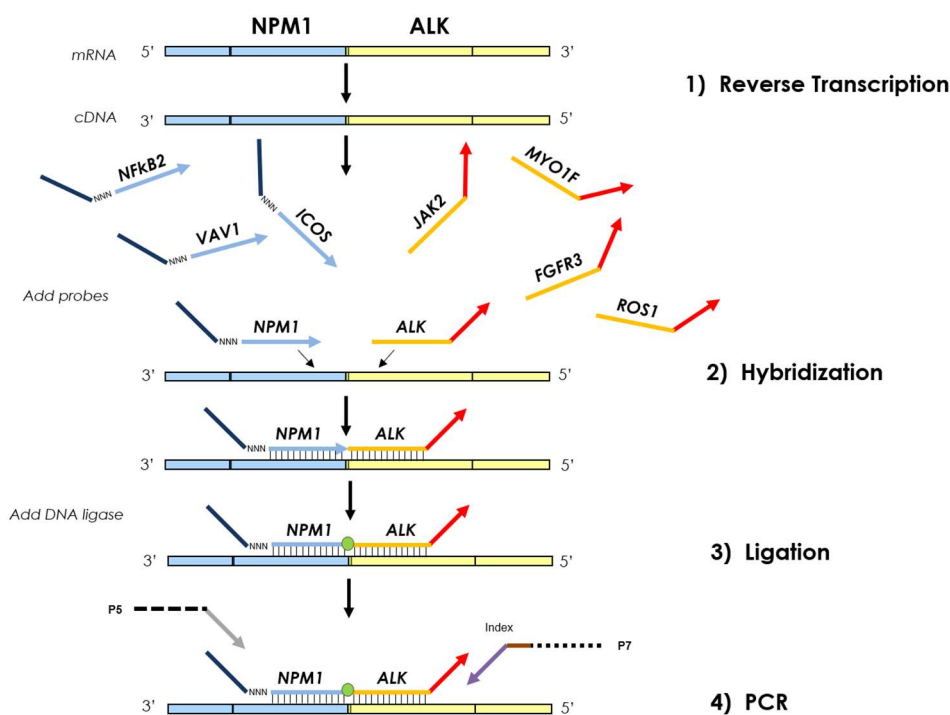
The fastQ files generated using this test contain data about counting sequences corresponding to the potential presence of a fusion transcript, i.e. ligation of two probes and their amplification.

They can be analysed using **GENEXPATH RT-MIS** software, which contains a specific sequence demultiplexing application.

By studying 69 genes, this test can detect fusion transcripts found in 7 types of T-cell non Hodgkin's lymphomas.


Test principle

The **GENEXPATH LymphoTranscript** test uses a ligation-dependent RT-PCR method (LD-RT-PCR). This semi-quantitative technique helps detect chromosome translations using specific oligonucleotide probe pairs. A probe pair for a control gene (GAPDH) is included in the test probe mix allowing an internal control of your experiment.

Four steps are sufficient to obtain libraries from a total RNA extraction.

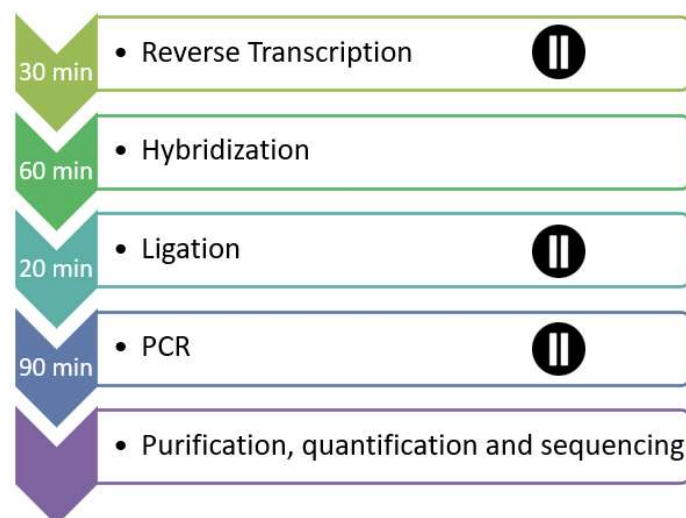- A reverse transcription (RT) step.
- A hybridization of specific oligonucleotide probes step.
- A ligation step.
- A PCR amplification step.

No purification is required until the libraries are obtained, which limits material losses and ensures this technique has excellent sensitivity. Also, the gene sequences targeted by the probes are particularly short (between 40 and 60 bases), which ensures excellent robustness with regards to RNA degradation.

LD-RT-PCR is therefore a particularly appropriate approach for analysing difficult biological samples like fixed paraffin-embedded tissue biopsies.

For each sample, around 10^5 sequences are sufficient to obtain an analysable expression profile, which helps test a large number of samples simultaneously in the same sequencing FlowCell. To optimise costs, **GENEXPATH LymphoTranscript** libraries can also be loaded at the same time as other sequencing libraries, generated by other methods.

Reagents

Content of GENEXPATH LymphoTranscript reagent kit

GENEXPATH LymphoTranscript probe mix	GEP-LT-PM
GENEXPATH LymphoTranscript barcodes	GEP-BC-xxx
GENEXPATH LymphoTranscript sequencing primer	GEP-SP-001
GENEXPATH LymphoTranscript GAPDH barcodes	GEP-BCC-xxx
GENEXPATH LymphoTranscript GAPDH sequencing primer	GEP-SP-002

XXX: barcode number

On receipt, these reagents should be stored between -30°C and -15°C.

They are ready to use and do not need to be diluted.

The shelf life of the reagents is 1 year.

Return to storage conditions immediately after use.

Do not use reagents after their expiration date stated on the label.

Format of reagent kits sold and quantities:

	Reagent kit - U = number of analyses		
	8U	16U	24U
GEP-LTPM probe mix	30 µL	48 µl	54 µl
Barcodes GEP-BC-xxx (from 033 to 044 depending on number of analyses purchased) BC=barcode	8 BC	8 BC	12 BC
	N°033 à 040	N°033 à 040	N°033 à 044
	5µL/BC	8µL/BC	6µL/BC
GEP-SP-001 sequencing primers	72 µL	120 µL	144 µL
For the internal control			
Barcodes GAPDH GEP-BCC-xxx (from 033 to 044 depending on number of analyses purchased) BCC=control barcode	8 BCC	8 BCC	12 BCC
	N°033 to 040	N°033 to 040	N°033 to 044
	5µL/BCC	8µL/BCC	6µL/BCC
GAPDH sequencing primers GEP-SP-002	72 µL	120 µL	144 µL

Reagents are supplied in larger quantities than actually required. After the ordered number of analyses has been completed, they should be discarded. If a new order is placed, new reagents will be delivered.

For a reagent kit with more than 8 analyses, each barcode will be used for 2 different analyses.

Reagents not supplied in the reagent kit:

Reagents	Suppliers and references
Invitrogen™ M-MLV Reverse Transcriptase (200U/µL)	Fisher Scientific, ref 10338842
Invitrogen™ dNTPs set (100mM)	Fisher Scientific, ref 10083252
Invitrogen™ Random Primers	Fisher Scientific, ref 10646313
SALSA MLPA Buffer (180 µL)	MRC Holland, ref SMR33
SALSA Ligase Buffer A (360 µL)	MRC Holland, ref SMR12
SALSA Ligase Buffer B (360 µL)	MRC Holland, ref SMR13
SALSA Ligase-65 (115 µL)	MRC Holland, ref SMR20
Q5® Hot Start High-Fidelity 2X Master Mix	New England BioLabs, ref M0494

AMPure XP (Magnetic beads)	Beckman Coulter, ref A63880
Qubit® dsDNA HS Assay Kit	Fisher Scientific, ref 10616763
Sequencing reagents	Illumina
TE buffer (10 mM Tris-Acetate pH 8.0, 1 mM EDTA)	Variable
Ethanol 100%	Variable
NaOH 1 N	Variable
Tris Buffer 200 mM pH 7	Variable
Nuclease-Free Water	Variable

On receipt and between each use, these reagents should be stored based on supplier recommendations.

Required equipment:

Equipment	Suppliers and references
Thermocycler in pre-PCR zone	Variable
Thermocycler in post-PCR zone	Variable
Qubit® Fluorometer (or equivalent)	Thermo Fisher Scientific, ref Q33238
Qubit® Assay Tubes	Fisher Scientific, ref 12037609
DynaMag™-96 Side Magnet - Magnetic plate (AMPure XP purification)	Thermo Fisher Scientific, ref 12331D
PCR 200 µL plates and tubes	Variable

Before starting

Biological samples

The **GENEXPATH LymphoTranscript** test is used to prepare sequencing libraries from total RNA extractions of tumour biopsies or human cell lines. This test is only applicable to Non-Hodgkin lymphomas.

These samples may be fresh, frozen or formalin-fixed, paraffin-embedded (FFPE).

To extract RNA from fixed tissues, we recommend using the kit Promega Maxwell® RSC RNA FFPE (Promega, ref AS1440 and AS4500).

The amount of RNA to be analysed should be between **50 and 500 ng, in a volume of 2 µL**. If the concentration of the solutions to be analysed is too high, this RNA can be diluted in nuclease-free water.

Solutions to prepare

- Random hexamers (Random Primers, Fisher Scientific, ref 10646313).
Working solution concentration: 100 µM
- Dilute 100 µL of the initial solution (at 3 µg/µL) by adding 1487 µL of nuclease-free water.
- Aliquot and store between -30°C and -15°C.
- dNTPs (dNTPs set, Fisher Scientific, ref 10083252).
Working solution concentration: 10 mM
- Mix the 4 stock solutions (250 µL each) and dilute by adding 1.5 ml of nuclease-free water.
- Aliquot and store between -30°C and -15°C.

Programming the thermocyclers

To limit the risks of contamination, use two thermocyclers, one in the pre-PCR zone and one in the post-PCR zone.

Two programmes are required:

- The first is for the first three steps of the protocol: **reverse transcription of RNA to cDNA, hybridization of oligonucleotide probes, and ligation**. It must be run on the thermocycler located in the pre-PCR zone.
- The second is used to **amplify ligation products and incorporate the barcodes and adapters required for sequencing**. It must be run on the thermocycler located in the post-PCR zone.

Programme 1: Pre-PCR

As the reaction volumes are small, ensure that the temperature of the thermocycler's heated lid remains high (95°C) at all steps of the programme to avoid evaporation.

Breaks at 4°C are provided between the different programme steps to add the required reagents.

Step 1: Reverse transcription.

Step 1a: Denaturation of RNA and hybridization of hexamers.

- Heated lid: 95°C
- 2 Minutes 80°C
- 5 Minutes 37°C

- 4°C infinite

Step 1b: Reverse transcription of RNA to cDNA.

- Heated lid: 95°C
- 15 Minutes 37°C
- 2 Minutes 98°C
- 4°C infinite

Step 2: Hybridization of probes.

- Heated lid: 95°C
- 2 Minutes 95°C
- 60°C infinite (1 hour of hybridization)

Step 3: Ligation.

- Heated lid: 95°C
- 54°C infinite (distribution of ligation mix)
- 15 minutes 54°C
- 5 minutes 98°C
- 4°C infinite

Programme 2: PCR.

- Heated lid: 95°C
- 6 minutes 94°C
- 35 x (30 seconds 94°C; 30 seconds 58°C; 30 seconds 72°C)
- 4 minutes 72°C
- 4°C infinite

Detailed protocol

Step 1: Reverse transcription

This step must be completed in the pre-PCR zone.

Required reagents.

- dNTPs 10 mM, random hexamers 100 µM, M-MLV RT kit (Reverse transcription buffer 5 x, DTT 100 mM, M-MLV RT enzyme), total RNA extraction to test (25 to 250 ng/µL).

It is recommended to carry out the entire procedure in 200 µL PCR plates or tubes.

Step 1.a: Reverse transcription; denaturation of RNA and hybridization of random hexamers.

- Thaw the following reagents, then keep them on ice or in a cooling rack:
 - Reverse transcription buffer 5 x
 - DTT 100 mM
 - dNTPs 10 mM

- Hexamers 100 μ M
- Prepare a reverse transcription mix. For each sample, mix (for a total volume of 3.75 μ L per reaction):
 - Reverse transcription buffer 1.25 μ L
 - DTT 0.5 μ L
 - dNTPs 1 μ L
 - Hexamers 1 μ L
- Distribute this mix in 200 μ L PCR tubes (3.75 μ L per tube) kept on ice or in a cooling rack:
- Add 2 μ L of each of the total RNA solutions (50 to 500 ng) to the different tubes.
- Vortex, centrifuge briefly.
- Place the tubes in the thermocycler in the pre-PCR zone, and proceed to **step 1.a of the Pre-PCR programme** (denaturation and hybridization of hexamers).

Step 1.b: Reverse transcription; conversion of RNA to cDNA

- At the end of step 1a, when the thermocycler temperature has dropped to 4°C, briefly centrifuge the tubes, then place them on ice or in a cooling rack.
- Add 0.5 μ L of reverse transcriptase (M-MLV RT) to each tube.
- Centrifuge briefly.
- Place the tubes back in the thermocycler.
- Check the temperature of the heated lid (95°C).
- Proceed to **step 1.b of the pre-PCR programme** (reverse transcription of RNA to cDNA).

Then proceed directly to step 2, or keep the ligation products between -30°C and -15°C.

Step 2: Hybridization of probes

This step must be completed in the pre-PCR zone.

Required reagents.

- **GENEXPATH LympHoTranscript** probe mix (GEP-LTPM), SALSA MLPA Buffer.

Hybridization of probes.

- **At the end of step 1b**, when the thermocycler temperature has dropped to 4°C, remove the tubes, centrifuge them briefly, and place them on ice or in a cooling rack.
- Thaw the Salsa MLPA buffer and the **GENEXPATH LymphoTranscript** probe mix, then keep them on ice or in a cooling rack.
- Prepare a hybridization mix. For each sample, mix (for a total volume of 3 µL per reaction):
 - Salsa MLPA Buffer 1.5 µL
 - **GENEXPATH LymphoTranscript** probe mix 1.5 µL
- Vortex, centrifuge briefly.
- Add 3 µL of this mix to each cDNA tube.
- Centrifuge briefly.
- Place the tubes back in the thermocycler.
- Check the temperature of the heated lid (95°C).
- Proceed to **step 2 of the pre-PCR programme** (hybridization of probes).

Step 3: Ligation

This step must be completed in the pre-PCR zone.

Required reagents

SALSA Ligase Buffer A, SALSA Ligase Buffer B, SALSA Ligase 65, nuclease-free water.

Ligation

- 15 minutes before the end of step 2a, thaw SALSA Ligase Buffer A and SALSA Ligase Buffer B and keep them on ice or in a cooling rack.
- Place the Salsa Ligase 65 enzyme on ice or in a cooling rack.
- Prepare a ligation mix. For each sample, mix (for a total volume of 32 µL per reaction):
 - Nuclease-free water 25 µL
 - Salsa Ligase Buffer A 3 µL
 - Salsa Ligase Buffer B 3 µL
- Vortex, centrifuge briefly

○ Salsa Ligase 65

1 µL

- Vortex, centrifuge briefly.
- After 60 minutes of incubation, proceed to **step 3 of the pre-PCR programme (ligation)**.
- Lower the temperature of the heated block to 54°C.
- Add 32 µL of the ligation mix directly to each tube, without removing them from the heated block.
- After distributing the mix, proceed to the next step of the programme (15 minutes at 54°C, 5 minutes at 98°C).

At the end of this step, when the temperature of the PCR block drops to 4°C, immediately proceed to step 4 (PCR amplification) or freeze the ligation products (between -30°C and -15°C).

After this step, do not keep products at higher temperatures (e.g. 4°C or room temperature) to avoid non-specific ligations which could result from residual enzyme activity.

Step 4: Amplification and incorporation of barcodes and adapters

In this step, the ligation products are amplified by PCR thanks to additional tails at the end of the probes. These amplifications are carried out using pairs of primers supplied in the **GENEXPATH LymphoTranscript** barcode tubes (GEP-BC-xxx).

To allow analysis of several samples in the same FlowCell, the PCR 3' primer has a molecular barcode which will be recognised by the demultiplexing algorithm of the **GENEXPATH RT-MIS** platform.

To carry out the internal control with GAPDH probes, two different PCR are completed, so you must double the number of tubes. For a given sample, you must use the same barcode number GEP-BC-xxx and GEP-BCC-xxx for information analysis. So you must add, for each sample, in one tube the barcode GEP-BC-xxx and in the other the associated barcode GEP-BCC-xxx.

Required reagents

GENEXPATH LymphoTranscript barcodes (GEP-BC-xxx), Q5® Hot Start High-Fidelity 2X Master Mix, nuclease-free water.

Amplification

- Prepare an amplification mix in the pre-PCR zone. For each sample, mix (for a total volume of 18 µL per reaction):
 - Q5® High-Fidelity 2X Master Mix 12.5 µL
 - Nuclease-free water 5.5 µL
- Vortex, centrifuge briefly.
- Distribute 18 µL of this amplification mix in the different wells in a PCR plate.
- Add 5 µL of ligation products generated in step 3 to each of the wells.
- Add 2 µL of **GENEXPATH LymphoTranscript** barcode (GEP-BC-xxx or GEP-BCC-xxx depending on test).

Use different GEP-BC-xxx barcodes for each tested sample, but for the same sample, use the same number for GEP-BC-xxx and GEP-BCC-xxx.

- Place the plate in the thermocycler in the **post-PCR zone**.
- Start **programme 2** (PCR).

At the end of the programme, when the temperature of the thermocycler drops to 4°C, quickly proceed to step 5 (purification) or freeze the amplification products between -30°C and -15°C.

Do not keep these products at higher temperatures for a long period (e.g. 4°C in the thermocycler or at room temperature).

Step 5: Purification and assay of sequencing libraries

At the end of the amplification step, the sequencing libraries should be purified to eliminate PCR primers and unincorporated nucleotides. This purification uses AMPure XP magnetic beads. The libraries should be assayed via fluorimetry before loading in the sequencer.

Required reagents

Ethanol 100%, nuclease-free water, AMPure XP beads, TE buffer (10 mM Tris-Acetate pH 8.0, 1 mM EDTA), Qubit® dsDNA HS Assay.

Step 5.a: Purification of sequencing libraries

Ensure that the beads are completely re-suspended before use.

- Purify 25 µL of PCR products with 45 µL of AMPure XP beads (following manufacturer recommendations).

- Elute the purified PCR products in 50 µL of TE buffer.

After purification, the libraries can be stored between -30°C and -15°C before sequencing.

Step 5.b: Assay of sequencing libraries:

- Assay 10 µL of each sequencing library via fluorimetry (Qubit dsDNA HS Assay kit, as per supplier recommendations).

Step 6: Dilution, pooling and sequencing of libraries

After purification, the **GENEXPATH LymphoTranscript** libraries must be diluted, pooled and loaded in the sequencer.

For optimal results, a minimum of 10⁵ sequences should be read for each sample.

Unlike most classic sequencing libraries, the reading of molecular barcodes required for demultiplexing **GENEXPATH LymphoTranscript** sequences takes place during read1. These sequences are therefore not demultiplexed automatically by the sequencer, and will be saved in “Undetermined” FASTQ files. Demultiplexing is carried out using the specific algorithm provided on the **GENEXPATH RT-MIS** platform.

Required reagents

- **GENEXPATH LymphoTranscript** sequencing primer (GEP-SP-001), **GENEXPATH LymphoTranscript** control sequencing primers (GEP-SP-002) (if internal control completed), Illumina sequencing reagents.

Sequencing on an Illumina MiSeq sequencer

For detailed information about dilution and denaturation of libraries, preparation of the sequencing primer, the injection sheet and the start of sequencing, refer to the Illumina guide to the MiSeq system.

- **Step 6.a: Dilution and pooling of libraries**

- Dilute each **GENEXPATH LymphoTranscript** library at a concentration between 2 nM and 4 nM, considering an average amplified fragment size of 150 pb.
- Pool **GENEXPATH LymphoTranscript** libraries in the equivalent volume.
- If other libraries are sequenced on the same FlowCell, adjust the concentrations of different pools, then combine them to obtain the desired sequence numbers (minimum 10⁵ sequences for each **GENEXPATH LymphoTranscript** library).

Example: For a pool of 10 **GENEXPATH LymphoTranscript** libraries requiring 1 M sequences (10⁵ sequences for each library), sequenced with a pool of libraries B at the same

concentration and requiring 3 M sequences, pool 1 μL of the **GENEXPATH LymphoTranscript** libraries pool and 3 μL of the pool of libraries B.

- *Step 6.b: Denaturation and dilution of the libraries pool*

- Denature and dilute the final pool based on recommendations in the Illumina guide to the MiSeq system, to obtain a final load concentration of 8 to 10 pM.

- *Step 6.c: Preparation of the sequencing primer*

- If the **GENEXPATH LymphoTranscript** libraries pool is sequenced alone, dilute 3 μL of each **GENEXPATH LymphoTranscript** sequencing primer (GEP-SP-001 and GEP-SP-002, if internal control) in a final volume of 600 μL of HT1 buffer, then place this 600 μL in well 18 of the MiSeq reagent cartridge.
- If the **GENEXPATH LymphoTranscript** libraries pool is loaded with other libraries sequenced using Illumina sequencing primers, pipette the entire content of well 12 (around 600 μL), add 3 μL of each **GENEXPATH LymphoTranscript** sequencing primer (GEP-SP-001 and GEP-SP-002 if internal control), then place this mix in well 18 of the cartridge.

- *Step 6.d: Preparation of the injection sheet*

- If the **GENEXPATH LymphoTranscript** library is sequenced alone, create the injection sheet to generate FASTQ files with 120 cycles in read 1.
- If **GENEXPATH LymphoTranscript** libraries are combined with other sequencing libraries, generate the injection sheet using the usual parameters, without entering the **GENEXPATH LymphoTranscript** samples.
- Specify the use of custom during run setup (with Local Run Manager on the Create Run page. In manual run mode, on the Run Setup screen).

In all cases, ensure that read 1 is carried out with a minimum of 120 cycles and the use of custom sequencing primer is specified.

- In all cases, the **GENEXPATH LymphoTranscript** library sequences will not be demultiplexed by the sequencer, but will be saved in an “Undetermined” FastQ file, which will then be loaded on the **GENEXPATH RT-MIS** platform.

- *Step 6.e: Start of sequencing*

- Start sequencing by following the procedure described in the Illumina guide to the MiSeq system.

Sequencing on an Illumina NextSeq 500/550 platform

For detailed information about dilution and denaturation of libraries, preparation of the sequencing primer, the injection sheet and the start of sequencing, refer to the Illumina guide to the NextSeq system.

- ***Step 6.a: Dilution and pooling of libraries***

- Dilute each **GENEXPATH LymphoTranscript** library at a concentration between 0.5 nM and 4 nM, considering an average amplified fragment size of 150 pb.
- Pool **GENEXPATH LymphoTranscript** libraries in the equivalent volume.
- If other libraries are sequenced on the same FlowCell, adjust the concentrations of different pools, then combine them to obtain the desired sequence numbers (minimum 10^5 sequences for each **GENEXPATH LymphoTranscript** library).

Example: For a pool of 10 **GENEXPATH LymphoTranscript** libraries requiring 1 M sequences (10^5 sequences for each library), sequenced with a pool of libraries B at the same concentration and requiring 3 M sequences, pool 1 μ L of the **GENEXPATH LymphoTranscript** libraries pool and 3 μ L of the pool of libraries B.

- ***Step 6.b: Denaturation and dilution of the libraries pool***

- Denature and dilute the final pool based on recommendations in the Illumina guide to the NextSeq system, to obtain a final load concentration of 0.8 pM to 1 pM.

- ***Step 6.c: Preparation of the sequencing primer***

- If the **GENEXPATH LymphoTranscript** libraries pool is sequenced alone, dilute 6 μ L of each **GENEXPATH LymphoTranscript** sequencing primer (GEP-SP-001 and GEP-SP002, if internal control) in a final volume of 2000 μ L of HT1 buffer, then place this 2 mL in well 7 of the MiSeq reagent cartridge.
- If the **GENEXPATH LymphoTranscript** libraries pool is combined with other libraries sequenced using Illumina sequencing primers, pipette the entire content of well 20 (around 2 mL), add 6 μ L of each **GENEXPATH LymphoTranscript** sequencing primer (GEP-SP-001 and GEP-SP-002 if internal control), then place this mix in well 7 of the cartridge.

- ***Step 6.d: Preparation of the injection sheet***

- If the **GENEXPATH LymphoTranscript** library is sequenced alone, create the injection sheet to generate FASTQ files with 120 cycles in read 1.

- If **GENEXPATH LymphoTranscript** libraries are combined with other sequencing libraries, generate the injection sheet using the usual parameters, without entering the **GENEXPATH LymphoTranscript** samples.
- Specify the use of custom during run setup (with Local Run Manager on the Create Run page. In manual run mode, on the Run Setup screen).

In all cases, ensure that read 1 is carried out with a minimum of 120 cycles and the use of custom sequencing primer is specified.

- In all cases, the **GENEXPATH LymphoTranscript** library sequences will not be demultiplexed by the sequencer but will be saved in the four “Undetermined” FastQ file, which will then be loaded on the **GENEXPATH RT-MIS** platform.
 - *Step 6.e: Start of sequencing*
- Start sequencing by following the procedure described in the Illumina guide to the NextSeq system.

Step 7: Results analysis

The sequence files generated by the Illumina sequencing platform (MiSeq or NextSeq) in FastQ format must be analysed using the **GENEXPATH RT-MIS** software available via the customer space at the following address: <https://connect.genexpath.com/>.

To help download the FastQ file, it should not be decompressed (fastq.gz).

This software is a comprehensive bioinformatics solution which includes different data processing algorithms. It carries out demultiplexing to assign sequences to each sample. It then precisely identifies the gene expression markers and quantifies them.

The **GENEXPATH LymphoTranscript** test is based on quantification of qualitative markers characterising the presence or absence of chromosome translocations.

GENEXPATH RT-MIS generates concise and transparent reports, from implementation of sequencing reactions to automated analysis of sequencing results.

GENEXPATH RT-MIS requires sequencing files to be loaded in FASTQ format, as well as the list of barcodes used during testing.

GENEXPATH RT-MIS assesses the sequencing quality of each sample by quantifying the number of reads identified and the number of UMI (unique molecular identifiers) detected.

For each sample, **GENEXPATH RT-MIS** generates an analysis report indicating the presence or absence of a fusion transcript, the number of reads and UMI obtained, as well as a bibliography reference corresponding to the transcript (if a fusion has been detected). This data can be downloaded.

GENEXPATH RT-MIS includes a user guide directly accessible online to explain how to use the tool, to describe all generated results and explain the presentation of results.

The company **GENEXPATH** does not permanently store results generated by the software **GENEXPATH RT-MIS**. The data must be downloaded directly after each analysis and saved by the user in their document management system.

Limits of the procedure

- The LymphoTranscript test was developed based on data from the literature to detect the most frequent fusion transcripts in patients with sarcoma. It is intended for FFPE use or frozen samples, possibly obtained from needle biopsies.
- A low amount of RNA or a low quality sample may cause an uninterpretable result.
- Sequencing must be performed with Illumina technology sequencers (Miseq and NextSeq).

Bibliography

[Detection of gene fusion transcripts in peripheral T-cell lymphoma using a multiplexed targeted sequencing assay.](#) Drieux et al., J Mol Diagn. 2021

Symbol table

 manufacturer	 reagent name
 manufacturing date	 temperature limit
 shelf life	 see instruction manual
 batch code	 For research use only

Notes

GENEXPATH LymphoTranscript reagents are protected by intellectual property rights and cannot be modified, reproduced, sold or transferred without the manufacturer's permission.

Information in this document is likely to change.

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